



EQAX-TC1

Trap Cartridge for CO₂

Specifications




Details	Specifications	
	<i>volume</i> ¹	<i>capacity</i> ²
1770-911, EQAX-TC1 Trap Cartridge for CO ₂	7 mL	≥ 200 L air
1770-912, EQAX-TC1 Trap Cartridge for CO ₂ (2-pack)	2×7 mL	≥ 400 L air

¹⁾ Volume of color-indicating carbon dioxide adsorbent based on functionalized macroporous polystyrene-divinylbenzene particles.


²⁾ Capacity was determined at continuous supply of carbon dioxide enriched air and then recalculated assuming a carbon dioxide concentration of 400 ppm in air.

Installation and use


Open storage bag, remove protective caps and connect cartridge to eluent bottle.

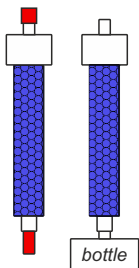

Attach protective caps when eluent is not withdrawn to maximize cartridge lifetime.

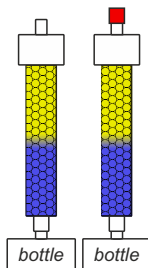

Stack cartridges when ~70% of capacity have been used to enable full capacity usage without risking eluent contamination.

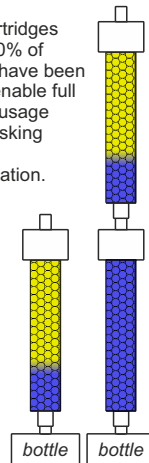

**Protect cartridges when not in use for longer lifetime.
Stack cartridges for full capacity usage.**

Xenoic[®] is a registered trademark of Diduco AB
Tvistevägen 48C, 90736 Umeå, Sweden
www.diduco.com · info@diduco.com